

# HIGH SENSITIVITY DOG CARDIAC TROPONIN-I ELISA

## Life Diagnostics, Inc., Cat. No. CTNI-4-HS

### INTRODUCTION

Cardiac troponin-I (CTNI) is a component of the troponin complex that regulates muscle contraction. After cardiac injury, CTNI is released into the blood. Because it is expressed specifically in the heart it is an excellent biomarker of cardiac injury. In humans, CTNI levels peak 12-24 hours after injury, returning to baseline within 2-6 days. In mice, levels peak as early as 1 hour and return to normal within 1-3 days.<sup>1</sup> Our ELISA kits have been used extensively since 2003 for measurement of CTNI in all species used in preclinical research. A representative list of publications citing their use can be found on our cardiac biomarker ELISA kits webpage (<https://lifediagnostics.com/cardiac-biomarker-elisa-kits/>).

### PRINCIPLE OF THE ASSAY

The ELISA uses two different antibodies that recognize a relatively protease-resistant epitope on CTNI. One is used for solid phase immobilization (microtiter wells). The second is conjugated to horse radish peroxidase (HRP) and used for detection. Standards and serum samples are incubated in the microtiter wells with HRP conjugate for one hour. This results in CTNI molecules being sandwiched between the immobilization and detection antibodies. The wells are then washed to remove unbound HRP-conjugate. TMB is added and incubated for 20 minutes. If CTNI is present a blue color develops. Color development is stopped by addition of Stop solution, changing the color to yellow, and absorbance is measured at 450 nm. The concentration of CTNI is proportional to absorbance and is derived from a standard curve.

### MATERIALS AND COMPONENTS

#### *Materials provided with the kit:*

- Anti-CTNI coated plate (12 x 8-well strips)
- CTNI Stock
- Diluent; YD12-1, 12 ml
- HRP Conjugate, 11 ml
- 20x Wash solution; TBS50-20, 50 ml
- TMB; TMB11-1, 11 ml
- Stop solution; SS11-1, 11 ml

#### *Materials required but not provided:*

- Pipettors and tips
- Distilled or deionized water
- Polypropylene or glass tubes
- Vortex mixer
- Absorbent paper or paper towels
- Plate incubator/shaker
- Plate washer
- Plate reader capable of measuring absorbance at 450 nm.
- Curve fitting software

### STORAGE CONDITIONS

Store the lyophilized stock at or below -20°C. The remainder of the kit should be stored at 2-8°C and the microtiter plate should be kept in a sealed bag with desiccant. Kits will remain stable for six months from the date of purchase.

### GENERAL INSTRUCTIONS

1. All reagents should be allowed to reach room temperature before use.
2. Reliable and reproducible results will be obtained when the assay is carried out with a complete understanding of the instructions and with adherence to good laboratory practice.
3. The wash procedure is critical. Insufficient washing will result in poor precision and falsely elevated absorbance readings.
4. Laboratory temperature will influence absorbance readings. Our ELISA kits are calibrated using shaking incubators set at 150 rpm and 25°C. Performance of the assay at lower temperatures will result in lower absorbance values.

### WASH SOLUTION PREPARATION

The wash solution is provided as a 20x stock. Prior to use, dilute the contents of the bottle (50 ml) with 950 ml of distilled or deionized water.

### STANDARD PREPARATION

1. Reconstitute the lyophilized CTNI stock with de-ionized or distilled water as detailed on the vial label. Mix gently until dissolved.
2. Label 7 polypropylene tubes as 10, 5, 2.5, 1.25, 0.625, 0.312, and 0.156 ng/ml.
3. Into the tube labeled 10 ng/ml, pipette the volume of diluent detailed on the stock vial label. Then add the indicated volume of stock and mix gently. This provides the 10 ng/ml standard.
4. Pipette 250 µl of diluent into the tubes labeled 5, 2.5, 1.25, 0.625, 0.312 and 0.156 ng/ml.
5. Prepare a 5 ng/ml standard by diluting and mixing 250 µl of the 10 ng/ml standard with 250 µl of diluent in the tube labeled 5 ng/ml. Similarly prepare the remaining standards by two-fold serial dilution.

The reconstituted CTNI stock should be frozen immediately after use. It remains stable when frozen or at least 1 month at -20°C and 6 months at -70°C. Discard the working standards after use.

### SAMPLE COLLECTION AND PREPARATION

Serum should be prepared as quickly as possible after blood collection and stored at 4°C. All samples should be similarly processed (i.e., storage times and temperatures should be the same). If serum samples cannot be assayed immediately they should be frozen at -70°C and thawed only once prior to use. Undiluted serum can be used with this kit. If dilution is necessary, use diluent YD12-1. Other diluents must not be used. Plasma cannot be used with this kit.

### ASSAY PROCEDURE

1. Secure the desired number of 8-well strips in the holder. Unused strips should be stored in the re-sealed bag with desiccant at 2-8°C for future use.
2. Dispense 100 µl of standards and samples into the wells (we recommend that standards and samples be run in duplicate).
3. Add 100 µl of HRP-conjugate into each well.
4. Incubate on a plate shaker at 150 rpm and 25°C for one hour.

5. Empty and wash the microtiter wells 5x with 1x wash solution using a plate washer (400  $\mu$ l/well).
6. Strike the wells sharply onto absorbent paper or paper towels to remove all residual droplets.
7. Dispense 100  $\mu$ l of TMB into each well.
8. Incubate on an orbital micro-plate shaker at 150 rpm and 25°C for 20 minutes.
9. After 20-minutes, stop the reaction by adding 100  $\mu$ l of Stop solution to each well.
10. Gently mix. It is important to make sure that all the blue color changes to yellow.
11. Read absorbance at 450 nm with a plate reader within 5 minutes.


### CALCULATION OF RESULTS

1. Using curve fitting software, construct a standard curve by plotting absorbance values of the standards versus  $\log_{10}$  of the concentration.
2. Fit the standard curve to a four-parameter logistic regression (4PL) equation (x axis =  $\log_{10}$  concentration) and determine the concentration of the samples.
3. Multiply the derived concentration by the dilution factor (if applicable) to determine the actual concentration in the original sample.
4. If the  $A_{450}$  values fall outside the standard curve, samples should be diluted appropriately and re-tested.

### TYPICAL STANDARD CURVE

A typical standard curve is shown below. This is for illustration only. A standard curve must be generated for each experiment.

CTNI (ng/ml)	$A_{450}$
10	2.736
5	1.362
2.5	0.796
1.25	0.448
0.625	0.280
0.313	0.194
0.156	0.138

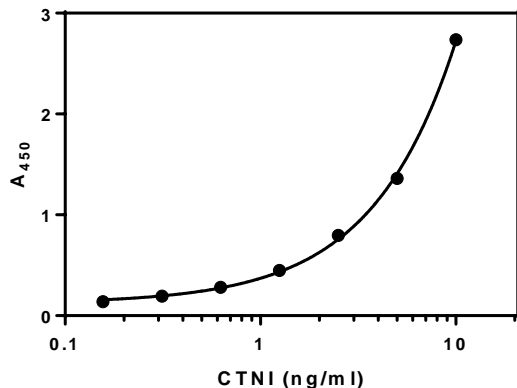

### REFERENCES

1. Engle SK. et al. Qualification of cardiac troponin I concentration in mouse serum using isoproterenol and implementation in pharmacology studies to accelerate drug development. Toxicologic Pathology. 37:617-628 (2009)

Rev 010518

For technical assistance please email us at  
techsupport@lifediagnosics.com