

Guinea Pig IgG ELISA

Life Diagnostics, Inc., Catalog Number: IGG-7

Guinea Pig IgG ELISA

INTRODUCTION

The guinea pig IgG ELISA kit is designed for measurement of IgG in guinea pig serum or plasma. The assay uses goat anti-guinea pig IgG for solid phase (microtiter wells) immobilization and horseradish peroxidase (HRP) conjugated goat anti-guinea pig IgG antibodies for detection. The immobilization and detection antibodies are cross-adsorbed with guinea pig IgA and IgM, ensuring specificity for IgG.

PRINCIPLE OF THE ASSAY

Test samples are diluted and incubated in the microtiter wells for 45 minutes alongside prepared guinea pig IgG standards. The microtiter wells are subsequently washed and HRP conjugate is added and incubated for 45 minutes. IgG molecules are thus sandwiched between the immobilization and detection antibodies. The wells are then washed to remove unbound HRP-labeled antibodies and TMB Reagent is added and incubated for 20 minutes at room temperature. This results in the development of a blue color. Color development is stopped by the addition of Stop Solution, changing the color to yellow, and optical density is measured spectrophotometrically at 450 nm. The concentration of IgG is proportional to the optical density of the test sample and is derived from a standard curve.

MATERIALS AND COMPONENTS

Materials provided with the kit:

- Anti guinea pig IgG coated 96-well plate (12 strips of 8 wells)
- Anti guinea pig IgG-HRP Conjugate, 11 ml
- Reference standard (lyophilized)¹
- 20x Wash Solution, 50 ml
- 10x Immunoglobulin Diluent, 25 ml
- TMB Reagent (One-Step), 11 ml
- Stop Solution (1N HCl), 11 ml

Materials required but not provided:

- Precision pipettes and tips
- Distilled or deionized water
- Polypropylene or glass tubes
- Vortex mixer
- Absorbent paper or paper towels
- Micro-Plate incubator/shaker mixing speed of ~150 rpm
- Plate washer
- Plate reader with an optical density range of 0-4 at 450nm
- Graph paper (PC graphing software is optional)

STORAGE OF TEST KIT

The unused kit should be stored at 2-8°C and the microtiter plate should be kept in a sealed bag with desiccant to minimize exposure to damp air. The test kit will remain stable for six months from the date of purchase.

¹ The reference standard consists of lyophilized guinea pig serum of known IgG concentration in a BSA matrix. The IgG concentration was determined relative to purified guinea pig IgG obtained from an independent laboratory.

GENERAL INSTRUCTIONS

1. Please read and understand the instructions thoroughly before using the kit.
2. All reagents should be allowed to reach room temperature (18-25°C) before use.
3. Optimum results are achieved if, at each step, reagents are pipetted into the wells of the microtiter plate within 5 minutes.

IMMUNOGLOBULIN DILUENT PREPARATION

The diluent is provided as a 10x stock. Prior to use estimate the final volume of diluent required for your assay and dilute one (1) volume of the 10x stock with nine (9) volumes of distilled or deionized water.

WASH SOLUTION PREPARATION

The wash solution is provided as a 20x stock. Prior to use dilute the contents of the bottle (50 ml) with 950 ml of distilled or deionized water.

STANDARD PREPARATION

1. The guinea pig IgG standard is provided as a lyophilized stock. Reconstitute as detailed on the vial label (***the reconstituted standard is stable at 4°C for one week but should be aliquoted and frozen at -20°C after reconstitution if future use is intended***).
2. Label 6 polypropylene or glass tubes as 100, 50, 25, 12.5, 6.25, and 3.13 ng/ml.
3. Into the tube labeled 100 ng/ml, pipette the volume of diluent detailed on the IgG standard vial label. Then add the volume of IgG standard indicated on the IgG standard vial label and mix gently. This provides the 100 ng/ml standard.
4. Dispense 250 µl of diluent into the tubes labeled 50, 25, 12.5, 6.25, and 3.13 ng/ml.
5. Prepare a 50 ng/ml standard by diluting and mixing 250 µl of the 100 ng/ml standard with 250 µl of diluent in the tube labeled 50 ng/ml.
6. Similarly prepare the 25, 12.5, 6.25, and 3.13 ng/ml standards by serial dilution.

SAMPLE PREPARATION

General Note: In our laboratory we have measured IgG levels in guinea pig serum of ~5 mg/ml. We suggest an initial dilution factor of 100,000 fold using the following procedure for each sample to be tested:

1. Dispense 498 µl and 798 µl of 1x diluent into separate tubes.
2. Pipette and mix 2 µl of the serum/plasma sample into the tube containing 498 µl of diluent. This provides a 250 fold diluted sample.
3. Mix 2 µl of the 250 fold diluted sample with the 798 µl of diluent in the second tube. This provides a 100,000 fold dilution of the sample.
4. Repeat this procedure for each sample to be tested.

Tissue extracts and tissue culture medium will likely have lower IgG levels than those found in serum. Optimal dilutions of such samples should be determined empirically.

ASSAY PROCEDURE

1. Secure the desired number of coated wells in the holder.
2. Dispense 100 μ l of standards and diluted samples into the wells (we recommend that samples be tested in duplicate).
3. Incubate on an orbital micro-plate shaker at 100-150 rpm at room temperature (18-25°C) for 45 minutes.
4. Aspirate the contents of the microtiter wells and wash the wells 5 times with 1x wash solution using a plate washer (400 μ l/well). The entire wash procedure should be performed as quickly as possible.
5. Strike the wells sharply onto absorbent paper or paper towels to remove all residual wash buffer.
6. Add 100 μ l of enzyme conjugate reagent into each well.
7. Incubate on an orbital micro-plate shaker at 100-150 rpm at room temperature (18-25°C) for 45 minutes.
8. Wash as detailed in 4 to 5 above.
9. Dispense 100 μ l of TMB Reagent into each well.
10. Gently mix on an orbital micro-plate shaker at 100-150 rpm at room temperature (18-25°C) for 20 minutes.
11. Stop the reaction by adding 100 μ l of Stop Solution to each well.
12. Gently mix. *It is important to make sure that all the blue color changes to yellow.*
13. Read the optical density at 450 nm with a microtiter plate reader within 5 minutes.


CALCULATION OF RESULTS

1. Calculate the average absorbance values (A_{450}) for each set of reference standards and samples.
2. Construct a standard curve by plotting the mean absorbance obtained from each reference standard against its concentration in ng/ml on linear graph paper, with absorbance values on the vertical or Y-axis and concentrations on the horizontal or X-axis.
3. Using the mean absorbance value for each sample, determine the corresponding concentration of IgG in ng/ml from the standard curve.
4. Multiply the derived concentration by the dilution factor to determine the actual concentration of IgG in the sample.
5. PC graphing software may be used for the above steps.
6. If the OD_{450} values of samples fall outside the standard curve, samples should be diluted appropriately and re-tested.

TYPICAL STANDARD CURVE

A typical standard curve with optical density readings at 450 nm on the Y-axis against IgG concentrations on the X-axis is shown below. This curve is for the purpose of illustration only and should not be used to calculate unknowns. Each user should obtain his or her data and standard curve in each experiment.

IgG (ng/ml)	Absorbance (450 nm)
100.00	2.7430
50.00	2.1050
25.00	1.3200
12.50	0.8050
6.25	0.4480
3.13	0.2740

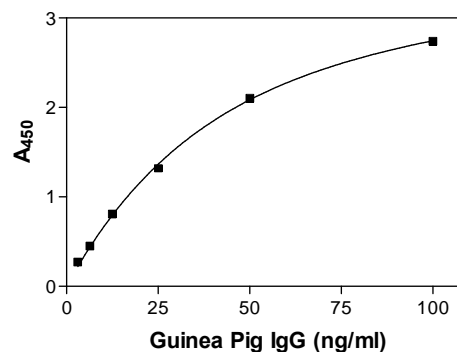

LIMITATIONS OF THE PROCEDURE

1. Reliable and reproducible results will be obtained when the assay procedure is carried out with a complete understanding of and in accordance with the instructions detailed above.
2. The wash procedure is critical. Insufficient washing will result in poor precision and falsely elevated absorbance readings.

Rev 041415NC

For technical assistance please email us at
techsupport@lifediagnosics.com