

USE STATEMENT

This kit is for research use only. Under no circumstances may it be used for diagnostic purposes.

INTRODUCTION

Attachment of polyethylene glycol (PEG) chains to therapeutic biologic agents, a process referred to as PEGylation, prolongs the circulating half-life of the modified protein by slowing proteolytic degradation and by masking it from the immune system. However, it has been reported that repeated injections of PEGylated proteins can induce anti-PEG antibodies that increase the rate of clearance and decrease drug efficacy (accelerated blood clearance, or ABC phenomenon). To aid research in this key area, we have developed a human anti-PEG IgA ELISA kit.

PRINCIPLE OF THE ASSAY

The assay uses immobilized mono-mPEGylated BSA (20 kDa PEG chain) as the capture antigen (coated on microtiter wells) and horseradish peroxidase (HRP) conjugated anti-human IgA heavy chain for detection. Serum or plasma samples are diluted and incubated alongside standards in the microtiter wells for 45 minutes. The wells are subsequently washed. HRP conjugate is added and incubated for 45 minutes. Anti-PEG IgA molecules are thus sandwiched between immobilized PEG and the detection antibody conjugate. The wells are then washed to remove unbound HRP-labeled antibodies. TMB reagent is added and incubated for 20 minutes at room temperature. This results in the development of a blue color. Color development is stopped by the addition of Stop Solution, changing the color to yellow. Optical density is measured spectrophotometrically at 450 nm. The concentration of anti-PEG IgA is proportional to the absorbance at 450 nm and is derived from a standard curve.

MATERIALS**Materials provided with the kit:**

- PEG-BSA coated plate (12 x 8-wells)
- Anti-IgA HRP Stock (lyophilized)
- Reference Stock (lyophilized)
- 20x HRP PEG Wash: PEGW50-20, 50 ml
- HRP PEG Diluent: PEGD50-1, 50 ml
- TMB: TMB11-1, 11 ml
- Stop Solution: SS11-1, 11 ml

Store at -20°C

Store at -20°C

Store at -20°C

Materials required but not provided:

- Pipettors and tips
- Distilled or deionized water
- Polypropylene or glass tubes
- Vortex mixer
- Absorbent paper or paper towels
- Plate incubator/shaker
- Plate washer
- Plate reader capable of measuring at 450 nm.
- Curve fitting software

STORAGE

The reference stock, HRP conjugate and the PEG-BSA coated plate should be stored at -20°C. All remaining kit components should be stored at 4°C. The microtiter plate should be kept in a sealed bag with desiccant. Kits will remain stable for six months from the date of purchase provided that the components are stored as described.

GENERAL INSTRUCTIONS/LIMITATIONS

1. Please read and understand the instructions thoroughly before using the kit.
2. All reagents should be allowed to reach room temperature (25°C) before use.
3. The wash procedure is critical. Insufficient washing will result in poor precision and falsely elevated absorbance readings.
4. Use only the wash solution and dilution buffer provided with the kit. PEG and PEGylated compounds are found in buffers conventionally used in ELISA's and cannot be used with this kit.
5. Kits are validated using plate shakers set at 150 rpm and 25°C. Performance of the assay at lower temperatures and/or mixing speeds will result in lower absorbance values.
6. Optimal results are achieved if at each step, reagents are pipetted into the wells of the microtiter plate within 5 minutes.

WASH SOLUTION

The wash solution is provided as a 20x stock. Prior to use, dilute the contents of the bottle (50 ml) with 950 ml of distilled or deionized water.

DILUENT

The diluent (PEGD50-1) is formulated for measurement PEG antibodies. It is supplied ready to use. DO NOT substitute other buffers.

STANDARDS

1. The anti-PEG IgA standard is provided as a lyophilized stock. Reconstitute as described on the vial label.
2. Label 8 polypropylene or glass tubes as 100, 50, 25, 12.5, 6.25, 3.13, 1.56 and 0 u/ml¹.
3. In the tube labeled 100 u/ml prepare the 100 u/ml standard as detailed on the stock vial label.
4. Dispense 250 µl of diluent into the remaining tubes.
5. Prepare a 50 u/ml standard by diluting and mixing 250 µl of the 100 u/ml standard with 250 µl of diluent in the tube labeled 50 u/ml.
6. Similarly prepare the 25 – 1.56 u/ml standards by serial dilution.

The reconstituted standard stock must be frozen at -20°C in a sealed vial if future use is intended.

SAMPLES

In studies at Life Diagnostics, Inc., we found that anti-PEG IgA levels in human samples ranged from undetectable to >23,000 u/ml. Optimal dilutions must be determined empirically. However, we suggest that samples initially be tested at 200-fold dilutions.

HRP CONJUGATE

Approximately 15 minutes before needed, reconstitute the lyophilized HRP conjugate as directed on the vial label and mix gently. Dilute as described on the vial label to give the working conjugate solution. The reconstituted conjugate stock must be frozen at -20°C in a sealed vial if future use is intended.

PROCEDURE

1. Secure the desired number of coated wells in the holder.
2. Dispense 100 µl of standards and diluted samples into the wells (we recommend testing in duplicate).
3. Incubate on a plate shaker at 150 rpm/25°C for 45 minutes.
4. Aspirate the contents of the microtiter wells and wash the wells five times with 1x wash solution using a plate washer (400 µl/well).
5. Strike the wells sharply onto absorbent paper to remove all residual wash solution.
6. Add 100 µl of diluted HRP conjugate into each well.
7. Incubate on a plate shaker at 150 rpm/25°C for 45-minutes.
8. Wash as detailed above.
9. Dispense 100 µl of TMB into each well.
10. Incubate on a plate shaker at 150 rpm/25°C for 20 minutes.
11. Stop the reaction by adding 100 µl of stop solution to each well.
12. Gently mix. It is important to make sure that all the blue color changes to yellow.
13. Read the optical density at 450 nm with a microtiter plate reader within five minutes².


RESULTS

1. Using curve fitting software, construct a standard curve by plotting absorbance values of the standards versus log₁₀ of the concentration.
2. Fit the standard curve using a second order polynomial (quadratic) equation and determine the concentration of the samples from the standard curve.
3. Multiply the derived concentration by the dilution factor to determine the actual concentration in the samples.
4. If the A₄₅₀ values of samples fall outside the standard curve, samples should be diluted appropriately and re-tested.

TYPICAL STANDARD CURVE

A typical standard curve with optical density readings at 450nm on the Y-axis against anti-PEG IgM concentrations on the X-axis is shown below. This curve is for the purpose of illustration only and should not be used to calculate unknowns.

Anti-PEG IgA (u/ml)	A ₄₅₀
100	3.258
50	1.742
25	0.936
12.5	0.513
6.25	0.268
3.13	0.172
1.56	0.118
0	0.067

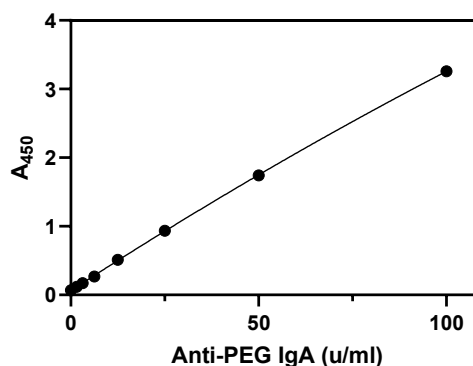


¹ One unit/ml is equivalent to 0.82 ng/ml of anti-PEG IgA


² On certain plate readers the A₄₅₀ value of the high standard may be out of range. If that occurs, absorbance values for all wells may be read at 405 nm instead. Absorbance values will be lower, but this does not affect results.

ASSAY PERFORMANCE

Parallelism: To assess performance of the assay, four samples sample containing anti-PEG IgA at concentrations of 725, 137, 62, and 3.5 u/ml were serially diluted to produce values within the dynamic range of the assay.

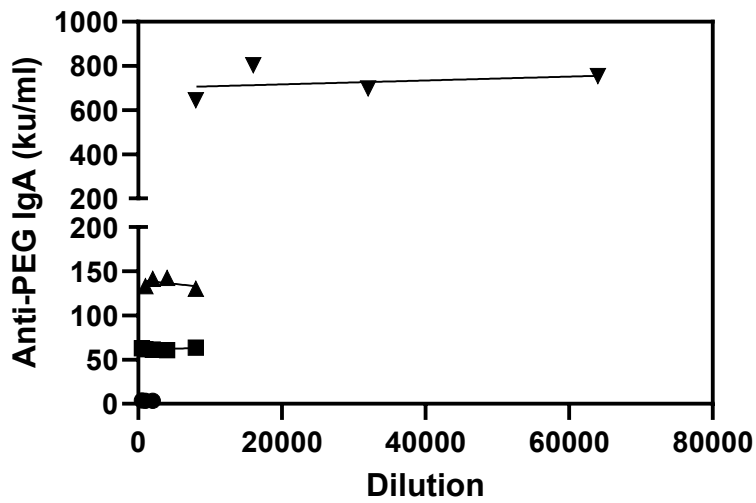

Anti-PEG IgA, IgG, and IgM: The heatmap shown below shows anti-PEG IgA, IgG, and IgM levels (ng/ml) in eighty human ascites fluid samples. The levels of anti-PEG IgA were higher than the levels of anti-PEG IgG or IgM.

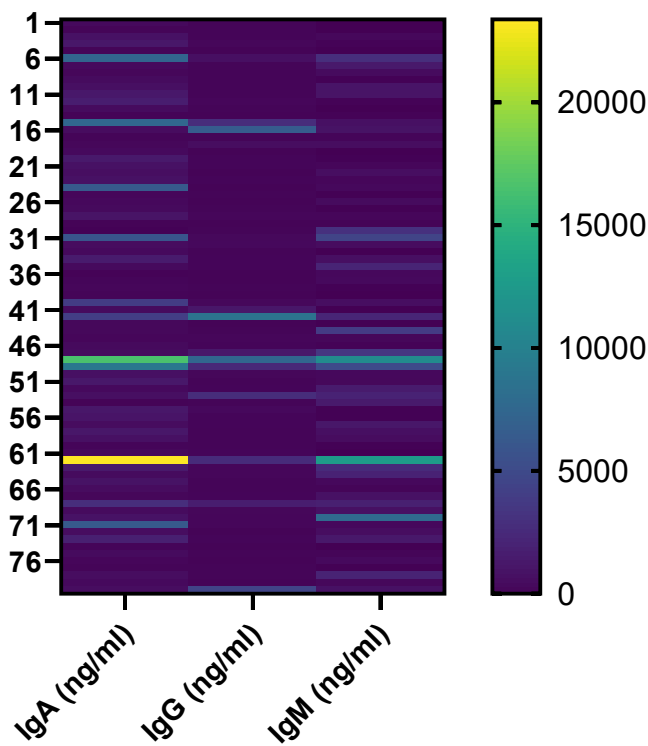

ASSAY UNITS

One unit/ml is equivalent to 0.82 ng/ml of anti-PEG IgA.

Rev 091625