

# COW SERUM AMYLOID A (SAA) SPARCL™ ASSAY

## Life Diagnostics, Inc., Catalog Number: SAA-SP-11

### INTRODUCTION

Serum amyloid A (SAA) is a positive acute phase protein of ~12 kDa that is expressed in the liver and circulates in blood. It is a useful biomarker of inflammation and disease in cows. Serum levels can increase >1500-fold during the acute phase response. In studies at Life Diagnostics, Inc., using this kit, we found SAA levels of  $11.1 \pm 13.1$  ng/ml in serum from healthy cows. Levels of  $2211 \pm 3415$  ng/ml and  $16,733 \pm 13,641$  ng/ml were found in serum from cows with mastitis and cows that were sick but non-mastitic, respectively<sup>1</sup>.

### PRINCIPLE OF THE ASSAY

The cow SAA SPARCL™<sup>2</sup> (Spatial Proximity Analyte Reagent Capture Luminescence, ref 1) assay uses two different peptide-specific cow SAA antibodies. One is conjugated to horseradish peroxidase (HRP), the other is conjugated to acridan, a chemiluminescent substrate. When the HRP and acridan conjugated antibodies bind to SAA they are brought into close proximity. With the addition of hydrogen peroxide, HRP catalyzes oxidation of proximal acridan molecules causing a flash of chemiluminescence. Acridan conjugated antibodies distant from HRP produce no signal. This principle allows the development of a homogeneous assay that allows rapid measurement of SAA concentrations.

The HRP and acridan conjugated antibodies provided with the kit are mixed with standards and diluted samples in wells of the 96-well SPARCL™ plate provided with the kit<sup>3</sup>. After incubation for 30 minutes on a shaker at 25°C and 150 rpm, the plate is placed into a luminometer. Trigger solution containing hydrogen peroxide is injected into each well and luminescence is immediately measured. The concentration of SAA is proportional to luminescence and is derived from a standard curve.

### MATERIALS AND COMPONENTS

- Anti-SAA HRP conjugate **Store ≤ -70°C**
- Anti-SAA acridan conjugate **Store ≤ -70°C**
- SAA stock **Store ≤ -70°C**
- Diluent; CSD50-1, 2 x 50 ml
- Trigger solution; TS7-1, 7 ml
- White SPARCL™ plate (12 x 8-well)
- Clear untreated 96-well plate

#### *Materials required but not provided:*

- Precision pipettes and tips
- Polypropylene microcentrifuge tubes
- Vortex mixer
- Plate incubator/shaker
- Luminometer capable of simultaneous injection/measurement
- Curve fitting software


### STORAGE

Store the HRP conjugate, acridan conjugate and SAA stock at -70°C (they may be stored at -20°C for one week). The remainder of the kit should be stored at 2-8°C. The SPARCL™ plate should be kept in a sealed bag with desiccant and antioxidant. The kit will remain stable

for at least six months from the date of purchase, provided that the components are stored as described.

### GENERAL INSTRUCTIONS

1. Please take the time to completely read all instructions before starting your assay. Contact us if you need clarification.
2. All reagents used in the assay should be allowed to reach room temperature (25°C) before use.
3. It is important that standards and samples be added to the SPARCL™ plate quickly. If testing large numbers of samples, rather than pipetting standards and samples directly into the white SPARCL™ plate using a single channel pipettor, we recommend the following. First, pipette an excess volume of standards and samples into appropriate wells of the clear 96-well plate. Then use an 8- or 12-channel multipipettor to quickly and efficiently transfer 50 µl aliquots to the appropriate wells of the white SPARCL™ plate. The wells of the clear plate have a maximum volume of 300 µl.
4. Follow the sequence of events below when running the assay.

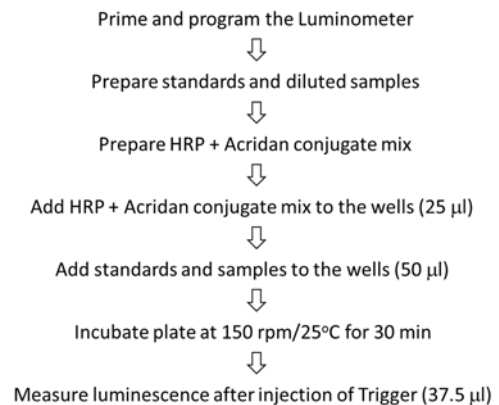

### STANDARD PREPARATION

The cow SAA stock is comprised of cow SAA lyophilized in a carrier protein matrix. The SAA content was determined by reference to a synthetic cow SAA polypeptide.

1. Reconstitute the stock with diluent as described on the vial label. Mix gently until dissolved.
2. Label 8 polypropylene tubes as 6.25, 3.125, 1.563, 0.781, 0.391, 0.195, 0.098 and 0.049 ng/ml.
3. Into the tube labeled 6.25 ng/ml, pipette the volume of diluent detailed on the stock vial label. Then add the indicated volume of stock and mix gently. This provides the 6.25 ng/ml standard.
4. Dispense 150 µl of diluent into the tubes labeled 3.125, 1.563, 0.781, 0.391, 0.195, 0.098 and 0.049 ng/ml.
5. Pipette 150 µl of the 6.25 ng/ml SAA standard into the tube labeled 3.125 ng/ml and mix. This provides the 3.125 ng/ml SAA standard.
6. Similarly prepare the remaining standards by two-fold serial dilution.

<sup>1</sup> The antibodies used in this kit are directed against the serum isoform of SAA, not the milk SAA isoform.

<sup>2</sup> The SPARCL technology was developed by Lumigen Corp.

<sup>3</sup> The plate provided with the kit has been treated with a reagent that reduces background chemiluminescence. Untreated plates cannot be used.

**Please Note: Use the standards within one hour of preparation.**

### SAMPLE PREPARATION

Serum or heparinized plasma should be prepared as quickly as possible after blood collection. If samples cannot be assayed immediately they should be frozen at or below  $-20^{\circ}\text{C}$ . Avoid repeated freeze-thaws.

The cow SAA SPARCL™ assay uses a homogeneous format and is therefore susceptible to a prozone or "hook effect" at high SAA concentrations. Samples should therefore be tested several dilutions in order to eliminate false low values. Because of the wide range of SAA concentrations that occur in serum throughout the acute phase response ( $\sim 10$  to  $>16,000$  ng/ml), testing over a range of dilutions also improves the chances of obtaining values within the standard curve range. We suggest testing dilutions of 20, 400 and 8000-fold:

1. Dispense 190  $\mu\text{l}$  of diluent into three separate tubes.
2. Pipette and mix 5.0  $\mu\text{l}$  of the serum sample into the first tube. This provides a 20-fold dilution.
3. Mix 10.0  $\mu\text{l}$  of the 20-fold diluted sample with 190  $\mu\text{l}$  of diluent in the second tube. This provides a 400-fold dilution.
4. Mix 10.0  $\mu\text{l}$  of the 400-fold diluted sample with 190  $\mu\text{l}$  of diluent in the third tube. This provides an 8000-fold dilution.

### CONJUGATE MIX PREPARATION

Instructions for preparation of the conjugate mix are detailed on the box that contains the HRP and acridan conjugates. If necessary, after thawing, briefly centrifuge to ensure that the contents are at the bottom of the tubes. Prepare the mix shortly before use using the diluent provided with the kit.

### LUMINOMETER SETUP

1. The luminometer must be capable of injection and simultaneous measurement of luminescence without any delay.
2. Prime the luminometer injection port with 1 ml of trigger solution.
3. Place the injection needle into the injection port as needed for BMG luminometers.
4. Program the luminometer to inject 37.5  $\mu\text{l}$  of trigger solution per well and to measure from time zero for 1 second (50 x 0.02 second intervals).
5. Define the format of the assay using the luminometer software.
6. Because the white SPARCL™ plate is provided as a 12 x 8-well strips, allowing use of fewer than 96-wells, make sure that the luminometer is programmed to inject trigger solution only into the wells being used.
7. We use a BMG LUMIstar Omega set at a gain of 3600. Optimal gain should be determined by the end user.
8. There are a number of manufacturers of luminometers that are equipped to run a SPARCL™ assay. Please contact Life Diagnostics or Lumigen ([www.lumigen.com](http://www.lumigen.com)) to discuss your luminometer.

### PROCEDURE

1. Before starting the assay ensure that the luminometer is primed with trigger solution and that the injection needle is positioned in the injection port.
2. Secure the desired number of SPARCL™ 8-well strips in the holder. Immediately seal unused strips in the resealable bag with desiccant and antioxidant. Store unused strips at  $2-8^{\circ}\text{C}$ .
3. Aliquot 25.0  $\mu\text{l}$  of conjugate mix into each well.
4. Dispense 50.0  $\mu\text{l}$  of standards and diluted samples into the wells (we recommend that standards and samples be tested in duplicate).

5. Incubate on an orbital micro-plate shaker at 150 rpm  $25^{\circ}\text{C}$  for 30 minutes.
6. After the 30-minute incubation, place the plate in the luminometer and measure luminescence after injection of trigger solution (37.5  $\mu\text{l}$ ).
7. Remove the plate from the luminometer and discard the used strips. Keep the plate frame if future use is intended.


### CALCULATION OF RESULTS

1. Before calculating results, review the raw data. If artefacts (RLU spikes) are apparent immediately after injection of trigger solution, eliminate that portion of the luminescence profile from analysis for all wells. We routinely use the sum of RLU values from a 100-980 ms data collection window.
2. Determine the sum of RLU values within the data collection window for the standards and samples.
3. Using graphing software, construct a standard curve by plotting the sum of the RLU values for the standards versus the  $\log_{10}$  of SAA concentration and fit to a sigmoidal, 4PL model.
4. Derive the corresponding concentration of SAA in the samples from the standard curve (remember to derive the concentration from the antilog).
5. Multiply the derived concentration by the dilution factor to determine the concentration of SAA in the original sample.
6. If the sum of the RLU values of diluted samples fall outside the standard curve, samples should be appropriately diluted and re-tested.

### TYPICAL STANDARD CURVE

A typical standard curve is shown below. This curve is for illustration only and should not be used to calculate unknowns. A standard curve should be run with each experiment.

SAA (ng/ml)	RLU
6.25	75580
3.125	38664
1.563	19090
0.781	6526
0.391	3785
0.195	2640
0.098	2005
0.049	1690

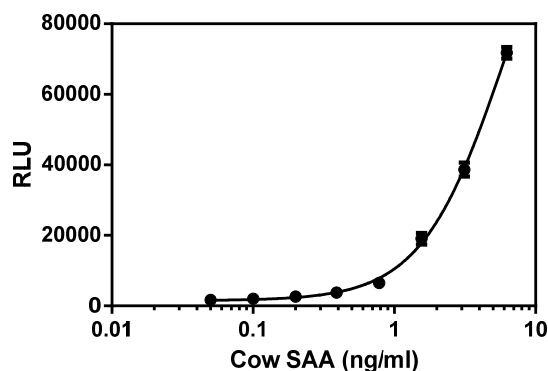

### REFERENCES

1. Akhavan-Tafti H. et al. A homogeneous chemiluminescent immunoassay method. *J Am Chem Soc.* 20;135(11):4191-4 (2013)

Rev 022118  
For technical assistance please email us at  
techsupport@lifediagnostics.com